



The FT90 Homogenisation Sub-System



The FT91 Homogenisation Sub-System

The FT90 and FT91 are complete in-line homogenisation sub-systems for use with Armfield miniature scale food processing equipment. The FT90 is based on an APV 1000 homogeniser and the FT91 is based on a Niro Twin Panda. Both homogenisers have been specifically optimised for this application. The units include controls, comprehensive instrumentation, valves, pipework, pulsation dampers, datalogging and all facilities required for operation and process integration with other Armfield equipment.

features

- Complete Homogenisation Subsystems
- Configurable for upstream or downstream homogenisation
- Variable throughput
- Two Stage Homogenisation with adjustable pressure (1000bar FT90, 600 Bar FT91)
- Twin Pistons (FT91) or Single Piston (FT90)
- Pulsation Dampers included
- Comprehensive instrumentation includes product temperature, product line pressure, homogenisation pressure and pump speed.
- All these parameters are displayed on the control console and available for datalogging via the integral USB datalogger
- Datalogging software seamlessly integrates the data with the data from the FT74 or similar.
- Adjustable divert facility for easy integration into continuous flow processes.
- Complete with interconnection hoses
- Can be sterilised and used in an Aseptic system

benefits

- In-line homogenisation for trial quantities and miniature-scale processing
- Simple integration with existing Armfield plant
- Uses scale-down valve technology which ensures the results are representative of production scale processing

Miniature-scale research & development technology





An ISO 9001 Company



Miniature-scale R & D technology

Head Office:

Armfield Limited
Bridge House, West Street,
Ringwood, Hampshire.
BH24 1DY England

Telephone: +44 1425 478781

Fax: +44 1425 470916

E-mail: sales@armfield.co.uk

U.S. Office:

Armfield Inc.
436 West Commodore Blvd (#2)
Jackson, NJ 08527
Telephone: (732) 928 3332
Fax: (732) 928 3542
E-mail: info@armfieldinc.com

Armfield R&D products include:

Hygienic connections
as standard



316 Stainless steel



CE certification



Extended warranty
as standard



For more product
information please visit:
www.explorearmfield.com



* Scan with smartphone with
QR code scanning software
installed.



* Scan for website



All products CE certified



follow us



twitter.com/Armfield_IFT
facebook.com/Armfielduk
linkedin.com/companies/armfield-limited
youtube.com/user/armfieldUK
explorearmfield.wordpress.com

0211/3k/S02796

Description

These units have been specifically developed to provide a simple and complete solution for in-line homogenisation in conjunction with the Armfield FT74 range of miniature scale processing equipment. They contain the homogeniser, pulsation dampers, a stainless steel control cabinet plus the required control valves, pipework and instrumentation all mounted within a stainless steel frame.

Hand wheel adjustment is provided for both stages of homogenisation pressure. The homogenisation pump speed is electronically controlled from the control console, and a manually adjustable divert valve is used to compensate for any inequalities between the flow rate from the FT74X and the homogeniser, or any fluctuations in these flows.

The homogenising valves, and the seats and balls of the pump valves, are manufactured from a hard wearing ceramic material, stellite.

The pulsation dampers are located upstream and downstream of the homogeniser to provide smooth flow on both sides. The pulsation dampers are non-pressurised and can be easily dismantled for cleaning. The stainless steel control console includes displays for the homogenisation pressure (both stages on FT90), for product temperature and for product process pressure. The outputs from all the sensors, together with the pump speed are also made available on an integral USB interface. The software provided with the FT90 and FT91 allows this data to be fully integrated with that from an FT74 datalogger, with a single display screen, single data table, and options to graphically display any or all of the data from either system.

The FT90 and FT91 are supplied with all necessary equipment, interconnecting hoses, homogeniser tools etc. for trouble free integration and operation.

Requirements - Services

Electrical Supply: (see Ordering Codes)
Water or Steam: for lubrication (FT90 only)
Steam: (sterile operation only)

Requirements - Accessories

A PC running Windows XP or later with an unused USB interface is required for datalogging. (If being used with FT74-44 or equivalent, two USBs are required).

For use with FT74-30MkIII, a homogeniser interconnection plate must be used. FT74-31 is used for upstream homogenisation (homogenisation during heating) and FT74-32 is used for downstream homogenisation (homogenisation during cooling)

Ordering Codes

FT90-E	380-415V/3ph/50Hz	3.5 kW
FT90-F	220V/3ph/60Hz	3.5 kW
FT91-A	220-240V/1ph/50Hz	2kW
FT91-G	230V/1ph/60Hz	2 kW

Shipping specification

	Gross weight	Volume
FT90	238kg	1.8m ³
FT91	240kg	1.8m ³

Overall dimensions

	H	W	D
FT90	1.15m	1.0m	0.95m
FT91	1.05m	1.1m	0.90m

Unit Specifications	FT90	FT91
Homogeniser	APV 1000	Niro Twin Panda 600
Throughput: Adjustable	8-22 L/hr	10-30 L/hr
Max 1st stage Homogenisation Pressures	1000 bar	600 Bar
Homogenisation Temperature	105°C max	90°C max
Max Sterilisation Temperature (30 mins)	130°C	140°C
Pump Power	3 kW	1.5 kW
No of pistons	1 off	2 off

The Armfield range includes HTST/UHT/aerobic systems, carbonator/filler/cappers, spray dryers/chillers, multifunction batch processors, ice cream freezers, margarine crystallisers, extractors, edible oils processors and more. For further information about our products and services, or to book a trial at one of our trials facilities, please contact us.

www.explorearmfield.com